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# EN - Safety cabinet operating instructions

## Chemicals, environmental, acid, alkali and gas cylinder cabinets

### 1. Intended application

All these cabinets are unsuitable for flammable liquids  
Cabinets for chemicals for storage of chemicals  
Environmental cabinets for storage of pollutant liquids  
Safety box - storing flammable liquids, aerosol cans etc.  
Cabinets for acids and alkalis - storage of aggressive liquids  
Gas cylinder cabinets - outdoor installation of gas cylinders

### 2. General safety notes

- Follow national/international laws and regulations for storage of hazardous materials and erection of safety cabinets
- Only use cabinet after obtaining proper instructions
- Any hazardous materials that escape must immediately be collected and removed.
- Follow limits (storage quantity, load etc.)
- Constructional changes to the product are not permitted

### 3. Erection / conditions for erection / functional checks

#### **All cabinet models**

Align cabinet squarely at its final location. Doors have to close and lock perfectly.

#### **Cabinets for acids and alkalis:**

- Attach anti-tilt bracket on the wall behind the cabinet. Place cabinet below the bracket
- If it is not possible to fasten to the wall, the cabinet must be fastened to the building using dowels through its floor.
- High cabinet with 4 adjusters (cabinet base). Adjust a max. of 3 vertically to a max. of 10 mm

#### **Gas cylinder cabinets:**

- Fix tightly to floor (2 holes,  $\varnothing$ 15 mm, base part) or wall (2 holes  $\varnothing$ 15 mm, upper rear panel)
- 4 adjusters(cabinet base). Adjust a max. of 3 vertically to a max. of 18 mm

### 4. Ventilation

Connection to a technical extraction air system is possible

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