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EN - Safety cabinet operating instructions

Chemicals, environmental, acid, alkali and gas cylinder cabinets

1. Intended application

All these cabinets are unsuitable for flammable liquids
Cabinets for chemicals for storage of chemicals
Environmental cabinets for storage of pollutant liquids
Safety box - storing flammable liquids, aerosol cans etc.
Cabinets for acids and alkalis - storage of aggressive liquids
Gas cylinder cabinets - outdoor installation of gas cylinders

2. General safety notes

- Follow national/international laws and regulations for storage of hazardous materials and erection of safety cabinets
- Only use cabinet after obtaining proper instructions
- Any hazardous materials that escape must immediately be collected and removed.
- Follow limits (storage quantity, load etc.)
- Constructional changes to the product are not permitted

3. Erection / conditions for erection / functional checks

All cabinet models

Align cabinet squarely at its final location. Doors have to close and lock perfectly.

Cabinets for acids and alkalis:

- Attach anti-tilt bracket on the wall behind the cabinet. Place cabinet below the bracket
- If it is not possible to fasten to the wall, the cabinet must be fastened to the building using dowels through its floor.
- High cabinet with 4 adjusters (cabinet base). Adjust a max. of 3 vertically to a max. of 10 mm

Gas cylinder cabinets:

- Fix tightly to floor (2 holes, \varnothing 15 mm, base part) or wall (2 holes \varnothing 15 mm, upper rear panel)
- 4 adjusters(cabinet base). Adjust a max. of 3 vertically to a max. of 18 mm

4. Ventilation

Connection to a technical extraction air system is possible

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